



**Networked Robotics Corp.**  
825 Chicago Ave, Suite F  
Evanston, IL 60202, USA  
Tollfree: (877) FRZ-TEMP  
(877) 379-8367

**General:** [info@networkedrobotics.com](mailto:info@networkedrobotics.com)  
**Technical:** [support@networkedrobotics.com](mailto:support@networkedrobotics.com)  
**Website:** [www.networkedrobotics.com](http://www.networkedrobotics.com)

---

## Release Notes NTMS Version 1.6 firmware

July 25, 2007

### ***About Firmware Upgrades***

Firmware upgrades are made to the NTMS by loading a Networked Robotics .nrf (Networked Robotics Firmware) file into the NTMS through the network using the Networked Robotics Hardware Upgrade Wizard. The Hardware Upgrade Wizard is available from the Download section of the Networked Robotics web site. Firmware upgrades can only be made from a PC on the same subnet as the NTMS.

If the network goes down, or a power failure occurs during the process of firmware upgrade, restart the NTMS unit and begin the process of firmware upgrade again.

### ***Compatibility***

NTMS firmware 1.6 can only be used with NTMS hardware shipped after November 1, 2006.

### ***Enhancements***

- 1) Support for the Dry Contact Probe (DCP)

This version supports a new Dry Contact Probe which can be used to connect the NTMS to devices with simple alarm contacts.